

# Technical Data Sheet

## Cured-Epoxy Remover - Formula 2

**#14942**

### Procedure

1. Dip slide with sections in solution A for 5 minutes.
2. Rinse slide in solution B - 2 changes of 1 minute each.
3. Rinse slide in solution C for 2 minutes.
4. Transfer slide to solution D for 1 minute.
5. Rinse in 100% ethanol - 2 changes of 1 minute each.
6. Rinse in 90% ethanol - 2 changes of 1 minute each
7. Rinse in 80% ethanol for 1 minute.
8. Bleach in solution E for 3 - 5 minutes.
9. Rinse again in 80% ethanol for 1 minute.
10. Wash with distilled water before staining.

Do not delay staining procedure once the resin has been removed from sections.